


Penlight Measuring Magnifier Battery and Bulb Replacement Instructions Product No. 753-1, 756

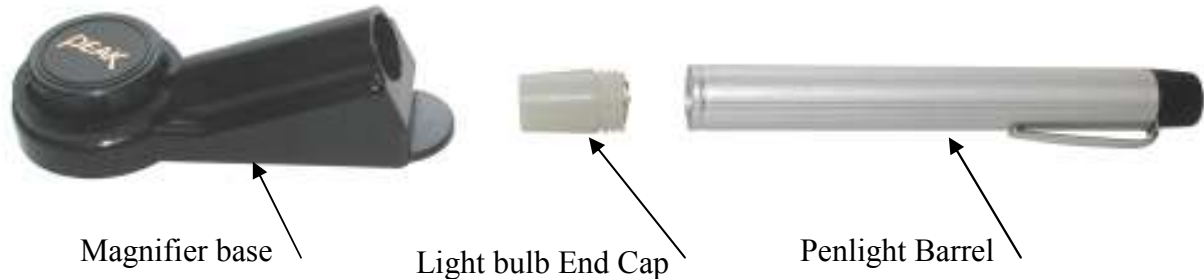

To install or replace batteries:

1. Pull the penlight barrel from the magnifier base. This is a press fit.
2. Unscrew the light end cap.
3. Insert two AA batteries with the positive terminal towards the light end cap.
4. Replace the light end cap.
5. Reinsert the penlight barrel in the magnifier until it seats on the bottom.
6. Turn on the light by twisting the rotary switch at the end of the penlight barrel.

To replace the penlight bulb:

1. Pull the penlight barrel from the magnifier base.
2. Unscrew the light end cap.
3. Remove the brass retaining ring and bulb
4. Replace lamp with a number 222 flashlight bulb (2.25V 0.25A).
5. Reinsert the brass retainer ring.
6. Insert two AA batteries with the positive terminal towards the light end cap.
7. Replace the light end cap.
8. Reinsert the penlight barrel in the magnifier until it seats on the bottom.
9. Turn on the light by twisting the rotary switch at the end of the penlight barrel.