

Form-ZERO™ (ZF)

Product Numbers 19986-10, 19986-11

Instructions for Use

For use with 10% Neutral Buffered Formalin (4% formaldehyde) and 4% Glutaraldehyde waste solutions. For use on aqueous solutions of 10% neutral buffered formalin and up to 4% glutaraldehyde.

1. Collect waste in the designated Waste Collection Container for Form-Zero™.
(1 gallon – Product Number 19986-13; 2.5 gallon – Product Number 19986-12)
2. Use a funnel with “screen insert” (Product Number 19986-7) while collecting waste to remove tissue debris from waste.
3. Add the entire contents of one Form-ZERO™ bottle (500ml) to each gallon (128-fl.oz.) or to 4-liters of aldehyde waste.
4. Secure lid on the Collection Container and agitate container to thoroughly mix solutions, making sure the agitation ultimately results in complete dissolution of Form-ZERO™.
5. Allow the mixed solutions to stand for 20 to 25 minutes for the reaction to be complete.
6. Before disposal, check with you company Safety Officer.
7. Pour treated waste into sanitary sewer and rinse container with cold tap water, preparing it for re-use.