

True North® Freezer Box Configurations

Product Numbers: 20962-1, 20962-2

INTRODUCTION

True North® Freezer Boxes are economical, reusable and moisture resistant plastic boxes for freezer storage. They are supplied as “flat-packed”, which eliminates the hassle of stacking and reduces the amount of space needed for storage. The clear plastic insert comes ready to hold tubes, cryovials, or mini centrifuge tubes (as shown in image on right); however, these inserts can be modified to also hold standard and large cryomolds. For examples of these configurations, see the sections below.



CONFIGURATION FOR STANDARD CRYOMOLDS



To configure the True North® Freezer Box for standard cryomolds, remove all vertical dividers from the insert except the 3rd and 6th one.

CONFIGURATION FOR LARGE CRYOMOLDS



To configure the True North® Freezer Box for large cryomolds, remove all vertical dividers from the insert except the 3rd and 6th one. Then, remove every other horizontal divider.

TED PELLA. INC.

Microscopy Products for Science and Industry

P.O. Box 492477, Redding, CA 96049-2477, U.S.A.

Telephone: 530-243-2200; 800-237-3526 (U.S.A. or Canada) • FAX: 530-243-3761

Email: sales@tedpella.com • Web Site: <http://www.tedpella.com>