

**STEREO VIEWER WITH CASE WITH 2X MAGNIFICATION  
PRODUCT NO. 758**



The #758 Stereo Viewer is a folding, pocket sized viewer which is ideal for three dimensional viewing of stereomicrographs. The precision ground and coated lenses give a two times magnification. A simple interpupillary adjustment calibrated in millimeters is provided. The viewer is packed in a protective carrying case.

To set up for viewing, simply pull down the folding legs. They will lock in place at the fixed focal distance. If needed, adjust the interpupillary distance. Place the stereo pairs side by side in the proper orientation for viewing. To do this, select the same spot or center area on each micrograph and slide and rotate (if necessary) gently until the stereo image appears. Keeping the micrographs flat greatly enhances the image.

Keep the stereo viewer clean and the lenses free from scratches by keeping it in its protective carrying case when not in use. Only use lens tissue for cleaning the lenses.