

FORELINE TRAP CONNECTION KIT FOR VRI-2 PUMP

92090-VRI

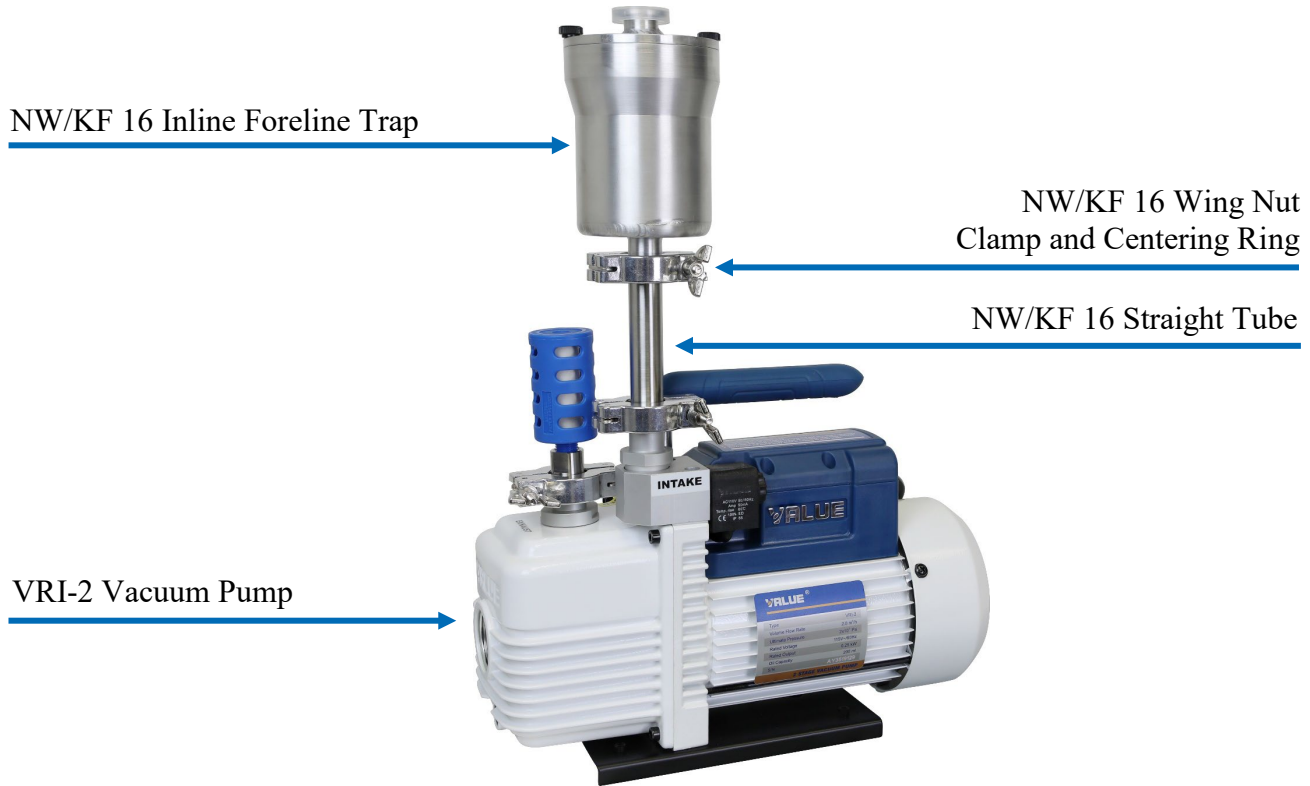


Figure 1: VRI-2 Rotary-Vane Pump with Foreline Trap (#92090-10) Installed

DESCRIPTION:

A vacuum connection kit to use the NW/KF 16 Inline Style Foreline Trap Products (#92090-10 and -30) with the VRI-2 Dual Stage Rotary Vacuum Pump Products (#92092 and 92092-230).

CONTAINS:

- 1x NW/KF 16 Straight Tube, 304 Stainless Steel (#92324-16)
- 1x NW/KF 16 Centering Ring, 304 SS with Viton® O-Ring (#92220-16)
- 1x NW/KF 16 Wing Nut Clamp, Cast Aluminum (#92210-16)

INSTALLATION:

1. Referencing **Figure 1**, use the Wing Nut Clamp and Centering Ring that were included with the VRI-2 Rotary-Vane Pump to attach the Straight Tube to the “INTAKE” on the pump.
2. Next, use the Wing Nut Clamp and Centering Ring included in the connection kit to attach the other end of the Straight Tube to the Inline Style Foreline Trap.